



## Starting/Stopping a WormBot experiment

1. Prepare the worms on a standard 12-well culture plate (this is covered in detail in a previous section/video)
2. Locate an empty slot(s) in the wormbot for your 12-well plate(s). The numbered plate locations for the WormBot are shown in the figure below:

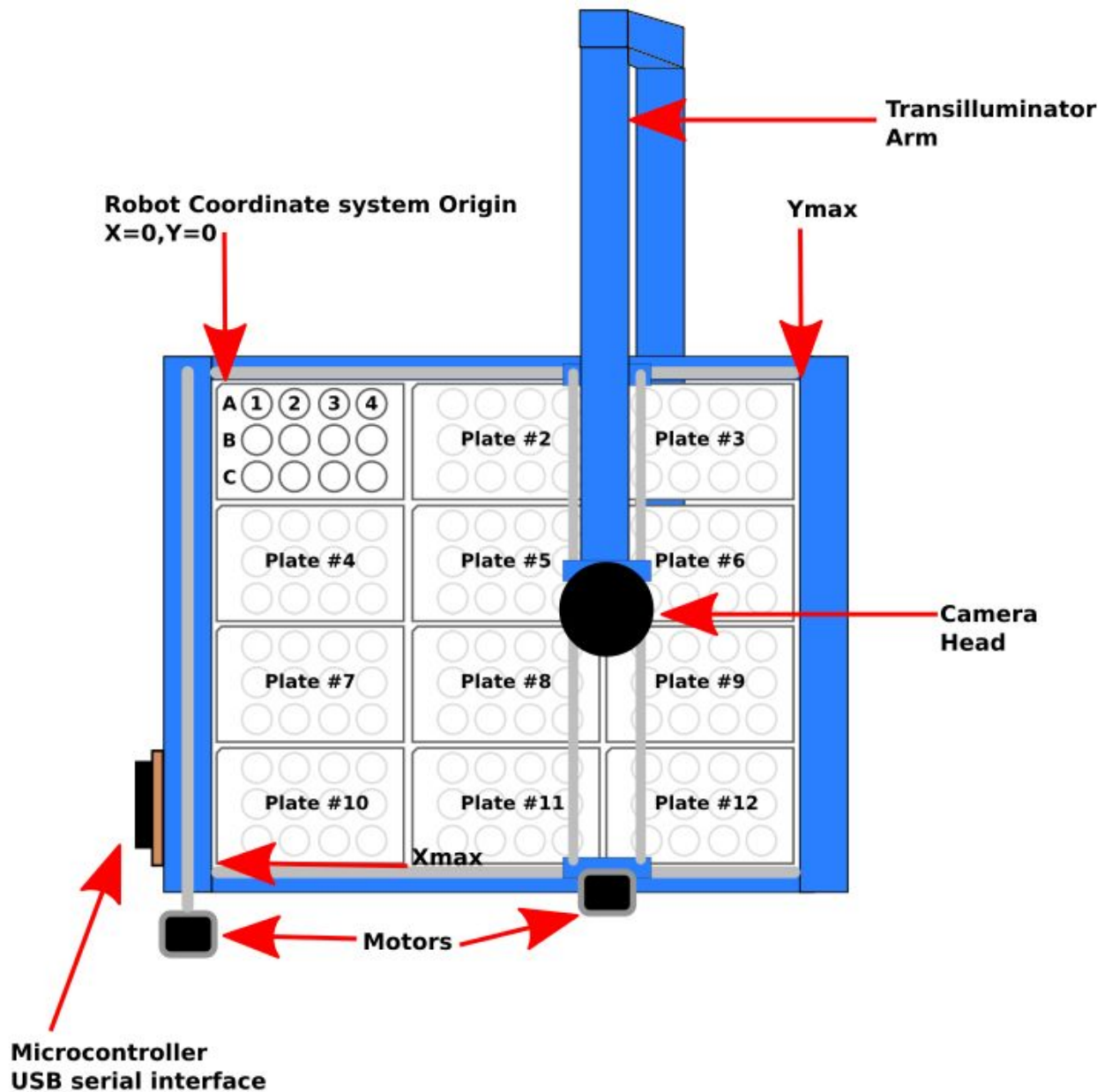

3. Open Firefox or another web browser on the server attached to your Wormbot and go to the Wormbot scheduler at <http://127.0.0.1/cgi-bin/scheduler> Note: You can get a static IP address or domain name for your WormBot so that you can access it from other locations. See the Security section for more details.

4. In the scheduler look at the list of available plate slots and select one from the drop down menu (labeled by red arrows below):

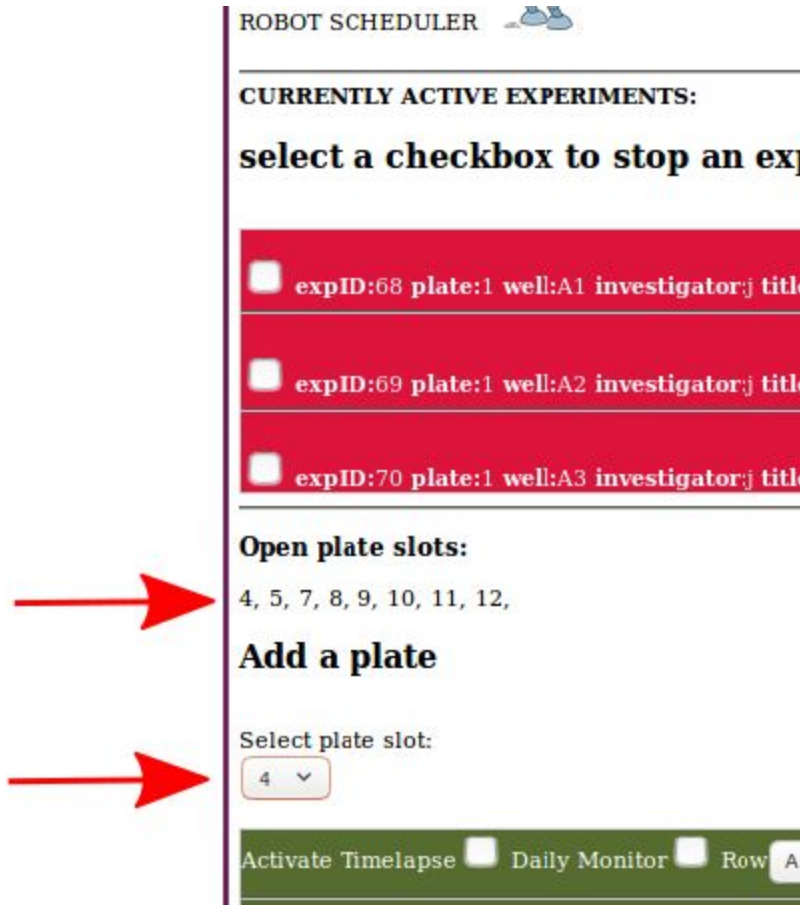

5. The green area of the scheduler is for inputting data about the experiments you are adding. There are 12 rows for the 12 wells in the plate (only one 12-well plate can be entered at a time). **Note: If you do not wish to collect data for all 12-wells you can leave some of the rows blank and data will not be collected.**

6. For each well there are 2 data collection modes available...both can be used simultaneously. If you select Activate Timelapse the WormBot will take an image of the well every 10 minutes until you stop the experiment. If you select Daily Monitor the WormBot will take a continuous video of the well (25 FPS) for 5 minutes once a day until you stop the experiment. Daily Monitor is useful for behavioral or other analyses that require frame rate faster than 1 frame every 10 minutes. While not necessary it is advisable to always Activate Timelapse data collection. Again, you may select both modes for a single well.

### Add a plate

Select plate slot:

Activate Timelapse <input checked="" type="checkbox"/>	Daily Monitor <input checked="" type="checkbox"/>	Row A
Activate Timelapse <input checked="" type="checkbox"/>	Daily Monitor <input checked="" type="checkbox"/>	Row A
Activate Timelapse <input checked="" type="checkbox"/>	Daily Monitor <input checked="" type="checkbox"/>	Row A


7. Enter the Row letter and well number for each Well/Experiment. **Note: the Green input window must be scrolled with the mouse wheel or window scroll bar to access all 12 wells.**

### Add a plate

Select plate slot:

Activate Timelapse <input checked="" type="checkbox"/>	Daily Monitor <input checked="" type="checkbox"/>	Row A	Well 3	Investiga
Activate Timelapse <input checked="" type="checkbox"/>	Daily Monitor <input checked="" type="checkbox"/>	Row A	Well 4	Investiga
Activate Timelapse <input checked="" type="checkbox"/>	Daily Monitor <input checked="" type="checkbox"/>	Row B	Well 1	Investiga

8. Enter the name of the investigator for each experiment and their email address and a title for the experiment and a description. As of the initial release the email field is not used however future releases of the Wormbot software will send experiment related emails to the address inputted here. **Note: certain ASCII characters are disallowed in these fields and you may notice them replaced in the final data files.**

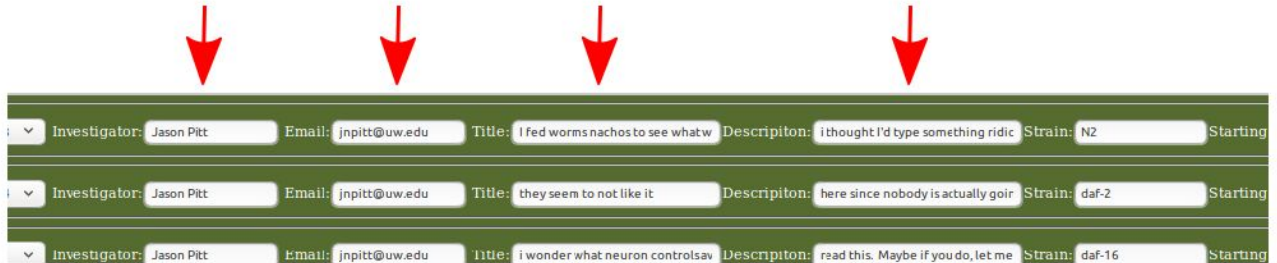

A screenshot of a data entry form with three rows. Each row contains fields for Investigator, Email, Title, Description, Strain, and Starting N. Four red arrows point downwards to the Title, Description, Strain, and Starting N fields of the first row.

Investigator:	Jason Pitt	Email:	jnpitt@uw.edu	Title:	I fed worms nachos to see whatw	Description:	i thought I'd type something ridic	Strain:	N2	Starting
Investigator:	Jason Pitt	Email:	jnpitt@uw.edu	Title:	they seem to not like it	Description:	here since nobody is actually goir	Strain:	daf-2	Starting
Investigator:	Jason Pitt	Email:	jnpitt@uw.edu	Title:	i wonder what neuron controlsav	Description:	read this. Maybe if you do, let me	Strain:	daf-16	Starting

9. Enter the starting N (number of worms placed into the well) and the age, in days, of the worms when the worms were placed on the Wormbot. **Note: As mentioned previously the best target N is somewhere between 30-40 animals and we usually transfer animals at L4 to the WormBot plates containing FUDR, however if you don't want to use FUDR you can transfer older worms after they've run out of sperm.**



A screenshot of a data entry form with three rows. Each row contains fields for Starting N and Starting Age(days). Two red arrows point downwards to the Starting N and Starting Age(days) fields of the first row.

Starting N:	33	Starting Age(days):	2
Starting N:	33	Starting Age(days):	2
Starting N:	33	Starting Age(days):	2


10. Once you've filled out all the fields, double check that all the data (especially the well letters and numbers) are correct press the Update Robot button.

### Add a plate

Select plate slot.

4 ▾

Activate Timelapse <input checked="" type="checkbox"/>	Daily Monitor <input checked="" type="checkbox"/>
Activate Timelapse <input checked="" type="checkbox"/>	Daily Monitor <input checked="" type="checkbox"/>
Activate Timelapse <input checked="" type="checkbox"/>	Daily Monitor <input checked="" type="checkbox"/>



The experiments should now show up in the Red Currently active Experiments section of the Scheduler:

---

**CURRENTLY ACTIVE EXPERIMENTS:**

**select a checkbox to stop an experiment**

<input type="checkbox"/>	expID:118 plate:4 well:A3 investigator:Jason Pitt title:I fed worms nachos to see what would happen i thought I'd type something ridiculous <a href="#">ANALYZE</a> <a href="#">VERIFY</a> <a href="#">STOP</a> <input type="checkbox"/>
<input type="checkbox"/>	expID:119 plate:4 well:A4 investigator:Jason Pitt title:they seem to not like it here since nobody is actually going <a href="#">ANALYZE</a> <a href="#">VERIFY</a> <a href="#">STOP</a> <input type="checkbox"/>
<input type="checkbox"/>	expID:120 plate:4 well:B1 investigator:Jason Pitt title:i wonder what neuron controls aver read this. Maybe if you doCOMMA let me know <a href="#">ANALYZE</a> <a href="#">VERIFY</a> <a href="#">STOP</a> <input type="checkbox"/>

---

**Open plate slots:**

11. To actually place the 12-well plate on the robot you can now simply put the plate into the empty slot you selected. However sometimes if the robot is busy this can be challenging or dangerous. After you have updated the robot, at the end of the next scan cycle (up to ten minutes) the robot will enter Loading Mode and the camera head will travel to the Xmax,Ymax position for a minute to give you time to load the WormBot while it is not scanning.

12. To stop an experiment simply check the two boxes per experiment line in the red Currently Active Experiments section of the scheduler. Then press the Update Robot button. As with starting an experiment, after the completion of the next scan cycle the robot will enter into Loading Mode, and the camera head will travel to Xmax,Ymax for a minute to allow you to unload the plate from the robot. **Note: You may stop individual wells on a plate and not others. Data will still be collected for experiments on that plate that have not been stopped. The plate slot will not become free to setup a new plate in the scheduler until all the wells in a plate have been stopped.**

**select a checkbox to stop an experiment**

<input checked="" type="checkbox"/>	expID:118 plate:4 well:A3 investigator:jason Pitt title:fed worras nachos to see what would happen i thought I'd type something ridi-ulous	<a href="#">ANALYZE</a>	<a href="#">VERIFY</a>	<a href="#">STOP</a>	<input checked="" type="checkbox"/>
<input checked="" type="checkbox"/>	expID:119 plate:4 well:A4 investigator:jason Pitt title:they seem to not like it here since nobody is actually going	<a href="#">ANALYZE</a>	<a href="#">VERIFY</a>	<a href="#">STOP</a>	<input checked="" type="checkbox"/>
<input checked="" type="checkbox"/>	expID:120 plate:4 well:B1 investigator:jason Pitt title:i wonder what neuron controls aver read this. Maybe if you doCOMMA let me know	<a href="#">ANALYZE</a>	<a href="#">VERIFY</a>	<a href="#">STOP</a>	<input checked="" type="checkbox"/>

**Open plate slots:**  
5, 7, 8, 9, 10, 11, 12.

**Add a plate**

Select plate slot:  
5

Activate Timelapse <input type="checkbox"/>	Daily Monitor <input type="checkbox"/>	Row A	Well 1	Investigator:	Email:	Title:	Description:
Activate Timelapse <input type="checkbox"/>	Daily Monitor <input type="checkbox"/>	Row A	Well 1	Investigator:	Email:	Title:	Description:
Activate Timelapse <input type="checkbox"/>	Daily Monitor <input type="checkbox"/>	Row A	Well 1	Investigator:	Email:	Title:	Description: